

# Validator<sup>®</sup> AVS Software 2.0

ACCURACY AND EFFICIENCY WITH SUPERIOR  
REPORTING AND STREAMLINED USER MANAGEMENT



In today's rapidly evolving digital landscape where Pharmaceutical and Biotechnology companies depend on accurate and reliable data, it is essential that reporting is streamlined and flexible. With the introduction of their latest software version 2.0 for Validator AVS, Kaye proudly presents a significantly enhanced reporting tool with simplified User Management capabilities.


The inclusion of Active Directory support in version 2.0, simplifies administrative tasks, enhances security,

and improves efficiency by providing a centralized platform for managing user authentication and access control.


Furthermore, the new software not only enhances software features, but it also enables connectivity to the renowned Kaye CTR-25 calibration bath. The CTR-25 liquid bath offers a volume of 2.5 liters, a temperature range of -25°C to 140°C, and has an exceptional stability of better than 0.01°C.

## FEATURES:


- Added Active Directory\* support
- Customized user types
- New user privileges added (Create, edit, and delete, split of Admin functions)
- New Common Reporting Tool integrated
  - Truncation of large studies to enable graph reports
  - Traceability between qualification studies and corresponding sensor calibrations
  - Pass/Fail report enhancements
  - Additional spread sheet report
- Additional audit trail report
- Parallel Report Generation with one click
- Generated Audit and Graph reports can be viewed in Reports Tile
- Improved data synchronization
- “Change console time” workflow modified to improve manual time sync
- Performance improvement during asset loading and navigating between different screens
- Support of the Calibration Bath CTR-25

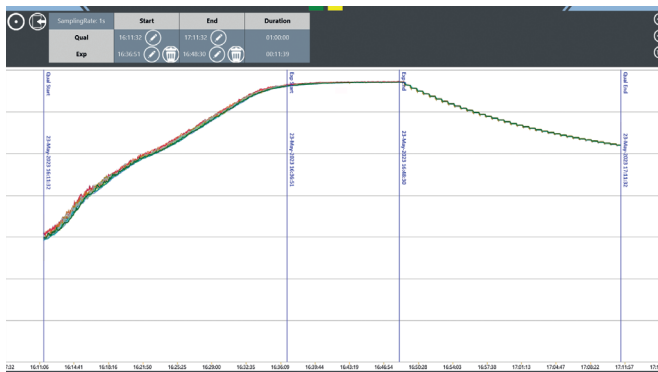


New Mark Cycles interface with enhanced workflow

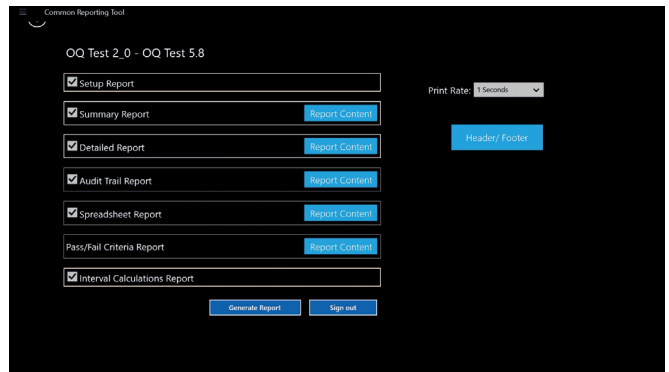

Generate all required reports with one click including audit trail data

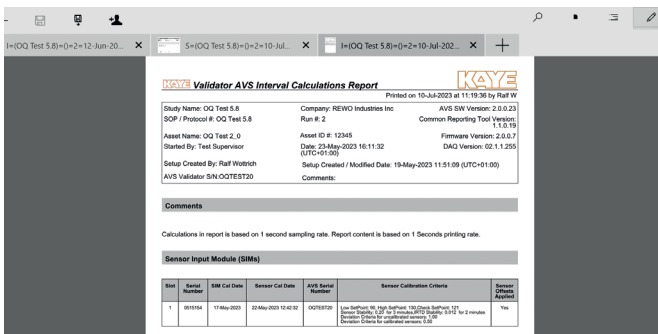

Enhanced traceability between sensor calibration

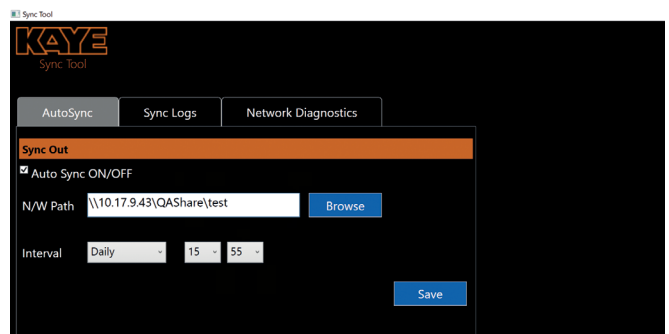

Sync Tool consistently running as a background service

## Kaye Representative contact:

**EUROPE, MIDDLE EAST, AFRICA AND ASIA**  
 Amphenol Advanced Sensors Germany GmbH  
 Sinsheimer Strasse 6  
 D-75179 Pforzheim  
**T:** +49 (0) 7231-14 335 0  
**F:** +49 (0) 7231-14335 29  
**Email:** kaye@amphenol-sensors.com

**USA/AMERICAS**  
 Amphenol Thermometrics, Inc.  
 967 Windfall Road  
 St. Marys, PA 15857  
**T:** +1(814) 834-9140  
**F:** +1(814) 781-7969  
**Email:** kaye-us@amphenol-sensors.com

## Request a demo:

**INDIA**  
 Amphenol Interconnect India Pvt Ltd.  
 Plot no. 6, Survey No.64 | Software Units layout  
 MAHAVEER TECHNO PARK  
 Hitech City, Madhapur | Hyderabad,  
 Telangana – 500081 | **T:** +91 40 33147100  
**Email:** kaye-india@amphenol-sensors.com

**CHINA**  
 Amphenol (Changzhou) Connector  
 Systems Co., Ltd, Building 10,  
 Jintong Industrial Park, No. 8 Xihu Road,  
 Wujin High-Tech Development Zone,  
 Changzhou, Jiangsu 213164  
**T:** 0086-519-83055197



**Warranty and disclaimer:** The information mentioned on documents are based on our current tests, knowledge and experience. Because of the effect of possible influences in an application of the product, they do not exempt the user from their own tests, checks and trials. A guarantee of certain properties or a guarantee for the proper suitability of the product for a specific, especially permanent application can not be derived from our data. Liability is therefore excluded to that extent permitted by law. Any proprietary rights of third parties as well as existing laws and regulations must be observed by the recipient of the product on his own responsibility.

© 2023 Amphenol Corporation. All Rights Reserved. Specifications are subject to change without notice. Other company names and product names used in this document are the registered trademarks or trademarks of their respective owners.